CF-10

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Table Top Centrifuge

Operating Manual

witeg Labortechnik GmbH

(English manual at page 10)

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Contents

1 Introduction	11
2 Safety Instructions	11
2.1 Description of symbols	11
2.2 Instructions.....	11
3 Product description	12
3.1 Key features.....	12
3.2 Control panel	13
4 Package content	13
5 Positioning the device and commissioning	13
6 Product Usage	14
6.1 Menu structure	14
6.2 Set Rotation Speed	14
6.3 Set delay time	14
6.4 SPIN DOWN	15
6.5 Key lock.....	15
6.6 Storage function.....	15
7 Instructions for Door Unlock	16
8 Troubleshooting Guide	16
9 Maintenance and cleaning	17
10 Technical data	17
11 Certificate	18

1 Introduction

Thank you for purchasing our **CF-10 Centrifuge**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a safe place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

**WARNING**

Information for avoiding injuries or fatal accidents.

**ATTENTION**

Safety instruction to avoid a damage of the product.

**ATTENTION**

Information about technical parameters.

**Environment**

Information about the operating environment.

**WARNING of explosions**

Safety instruction to avoid explosions.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.
- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.

- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.

- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- All circuits and electrotechnical components are **witeg** owned technology. In fact of this it is necessary that only **witeg** authorized personnel repair the device.
- Pull out the power cable if you don't want to use the device for a longer time.

- It is important that the device stands up on a horizontal, stable and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Repairs and amendments should only be made by **witeg** authorized people. Contraventions exclude the warranty.

- Do not use this device in areas with potentially explosive atmosphere.


3 Product description

3.1 Key features

- Only 57 dB at maximum speed.
- Maintenance-free Motor.
- High performance (ca. 12.200 x g).
- Compact and durable design.
- Simplest operation by patented Jog-Shuttle controller.
- Illuminated LCD with power bar graph.

- Acceleration and breaking time lower than 15 seconds.
- For 12 x 0,2-/0,5-/1,5-/2,0 ml tubes.

3.2 Control panel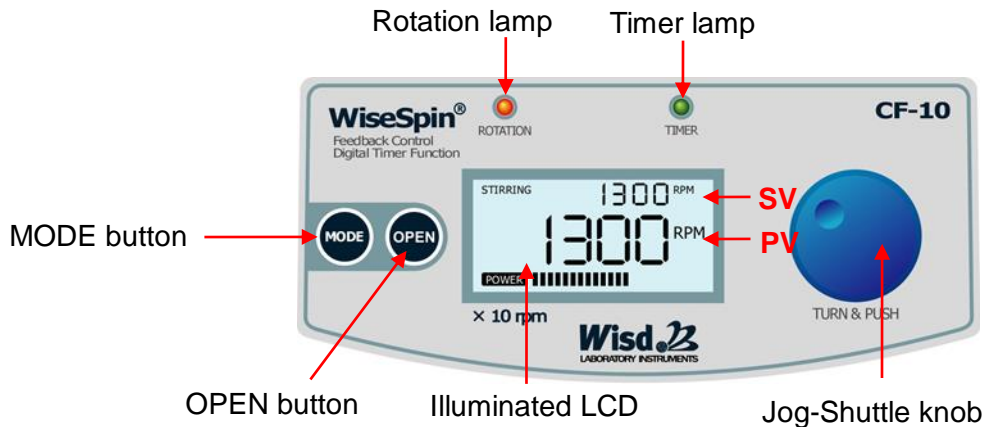

4 Package content


quantum	denomination
1	CF-10 Table Top Centrifuge
12	0.2-/0.5- ml PCR tube Adaptor
1	power supply curt
1	operating manual

5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol: 
2. Remove the package carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your costumer service.
4. Remove perhaps soiling and package material and other items in the centrifugal chamber.
5. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
6. Make sure that you follow all safety instructions.
7. Connect the device with the power outlet and Turn on the Main Switch.
8. The Air hole is closed with rubber plug for shipment at factory. Remove the plug for use.
9. Push the "Open" button to open the door and check the rotor for abnormality. And, put the sample tube of same weight for symmetry.
10. Press strongly the Door (Lid) so that it is completely closed with the Electronic Safety Lock.


6 Product Usage

6.1 Menu structure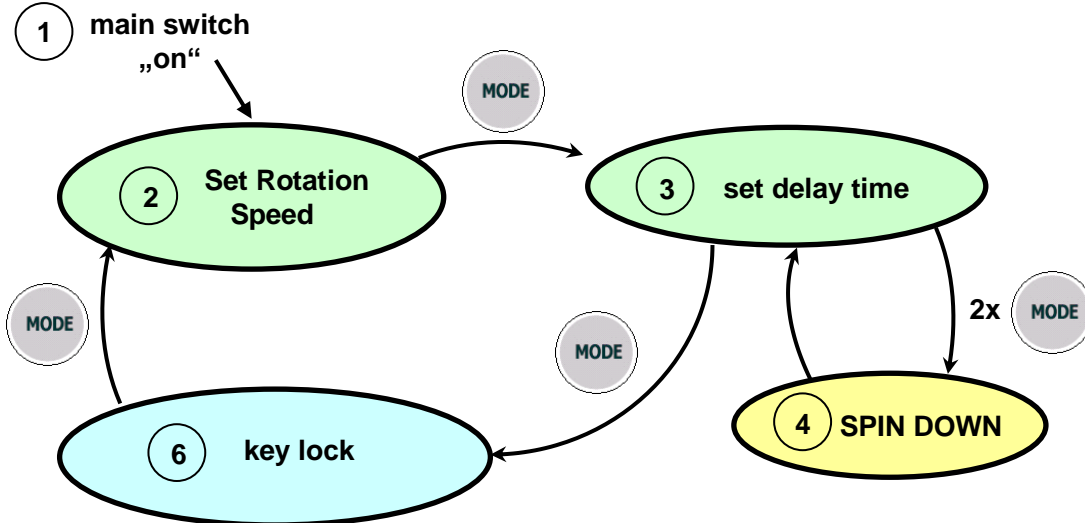

6.2 Set Rotation Speed


- When the power is turned ON, the initial screen is the setting mode of the set Rotation speed mode.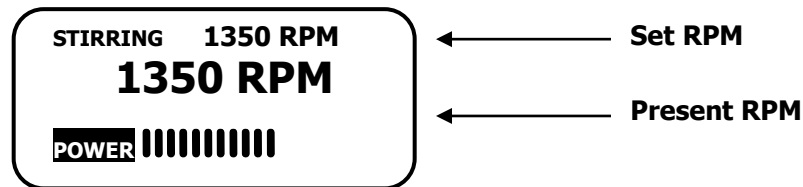

- ROTATION LED is lighted during the operation of Rotation, and indicates that it is operating.
- If “STIRRING” is blinking on the LCD there is a speed value stored. The saved value is showed in the SV you can start the stirring process by pushing the Jog-Shuttle knob.
- In both cases is the speed during the stirring process variable. To change the SV you only have to turn the Jog-Shuttle knob clockwise to increase the speed and counter-clockwise to reduce it.
- To stop the stirring process manually, you have to push the Jog-Shuttle knob again or decrease the speed till “0” by turning the Jog-Shuttle knob counter-clockwise.
- Although the Stop button and the Open button are pushed during operation, the door does not open until the Rotator stops completely. It is for safety. When the Rotator completely stops, the Door opens automatically.
- It takes only less than 15 seconds for the Rotator of the highest speed to stop completely. It also takes only less than 15 seconds for the stopped Rotator to reach the highest rotation speed.

6.3 Set delay time

- If pressing the MODE button in the Set Rotation Speed Mode, it is switched to the Set Timer mode.
- TIMER LED is lighted during the operation of Timer, and indicates that it is operating.
- Timer will automatically run, when after reach to set RPM at Set RPM mode.

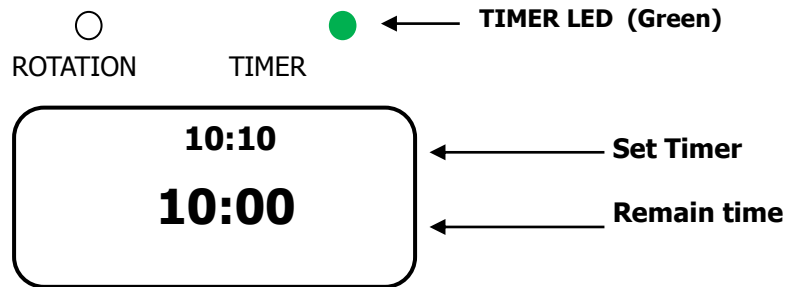- The range of setting time is from 00:00 to 99min and 59sec, and 00:00 indicates continuous operation.
- You can change or save the setting by using the operation method of the Jog-Schuttle, and Start/Stop of operation can be performed.

6.4 SPIN DOWN

- This mode is very convenient for repetitive use of certain rotation speed..

- In the timer mode, a further turn of the Jog-Schuttle Knob from the timer setting of "0" to the left indicates "SPIN" on the LCD screen to enable the Spin Down function.
- Push the Mode button twice to switch to the Set Rotation Speed mode, and then set the Spin Down RPM.
- When the Spin Down RPM starts to be activated, it reaches the set RPM and then immediately slows down and stops.
- When other timer function than the Spin Down is needed later, make sure to set the timer at higher than "0".
- You can change or save the setting by using the operation method of the Jog-Schuttle, and Start/Stop of operation can be performed.

6.5 Key lock

- To avoid accidental changing of the set values, push the MODE button in the menu item "set delay time".
- In this menu item can you see the actually speed and Time. The function of the Jog-Shuttle knob is deactivated.
- Press the MODE button once more to leaf the key lock. Than you are in the menu item "set speed".

6.6 Storage function

The **CF-10 Table Top Centrifuge** includes a storage function for speed and delay time values. That is practical if you want to work with the same parameters often. Then you need not to set them in every usage.

- For storing values you only need to hold the Jog-Shuttle knob pushed for two seconds. Successful storing is displayed by blinking of the SV for three times.
- If you saved a speed value, "STIRRING" is blinking and after the devices restart.
- If you saved the delay time it is not showed at the LCD.
- Every stored parameter is erasable. Turn the value with the Jug-Shuttle till "0" and then hold the knob pressed for two seconds.

- If you want to store a new value it is not necessary to erase the old one.

7 Instructions for Door Unlock

- When this equipment operates the centrifugal separation, its door is designed not to be opened even with the “OPEN” button for the purpose of safety. Door opens is only after centrifugation is completed. But the door can open manually in the event that it is to be inevitably opened while the power is off or it cannot be opened with the faulty “Open” button.

- As shown in the figure, the door can open manually by inserting a thin long tool such as clip tip into the hole of keyway.
- Try it only when the power is off or the open button is faulty. The door can open manually even while centrifugal separation is operated, but it is allowed only in unavoidable situation in which the operation cannot but stop forcibly for some reasons such as incorrect mounting of sample

* (NOTE) If you open the lid during rotation, the unit will automatically stop the operation for user safety. Do not open the door during rotation !! If you don't keep this note, you can have a serious problem of the machine, and so it leads to an accident as well.

8 Troubleshooting Guide

Situation	Confirmations and Solutions
Power Failure	<ul style="list-style-type: none"> ✓ Check the supplied voltage is present. ✓ Check the power connection and the power cord for proper connection and integrity. ✓ Check if the MAIN power switch is ON.
Not Rotation	<ul style="list-style-type: none"> ✓ Check the set RPM. ✓ Check the Lid is opened. ✓ Check the any prevent material beside rotator. ✓ Check if the MAIN power switch is ON.
ERROR 1	<ul style="list-style-type: none"> ✓ Check the set RPM and set timer.

If other problems arise, or one of the above not solves problems, please contact your official agent or the manufacturer.

Repairs and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

9 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts and accessories.

10 Technical data

Model	CF-10
Dimensions (mm) (w×d×h)	170 x 226 x 183
Speed (rpm)	1.000 – 13.500
Display	Illuminated LCD
Capacity	12 x 0,2-/0,5-/1,5-/2,0 ml tubes
Overall power consumption	90 W
Accelerating time	15 s to max. speed
Stop time	20 s from max. speed
Motor	Brushless DC motor, direct drive
Sound level	≤ 57 dB
Weight (kg)	2,3
Safety devices	Auto stop, Motor-error sensor
Power supply	AC 230 V, 1N~, 50/60 Hz

11 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	CF-10 Table Top Centrifuge
Serialno.	
Date	

witeg Labortechnik GmbH

Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de